


LAE 3 Fridge Controller Guide

This display shows the temperature. This reading is showing 3°.

When the light is showing below the snowflake, this means the compressor is running.

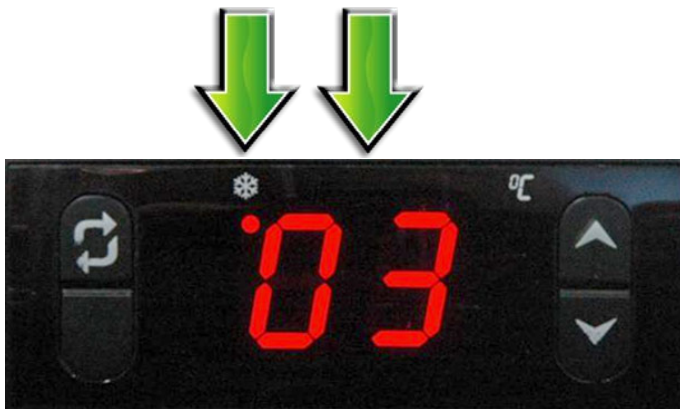

This button is to check the low temperature.



This low temperature is reading -1°

This button is to check the high temperature.




This high temperature is reading 12°

When the display reads “AL” this means the high or low temperature has been surpassed. To silence the alarm press the up/down buttons. The current temperature may also display between the AL flashing on the screen.



Press the up/down arrows to check why it alarmed before resetting the memory.

Resetting the memory will remove ALL the data stored in memory, including the high & low temperature and alarm readings. To reset the memory, **press and hold** the top left button with the two chasing arrows for 3 seconds until “rr” is displayed.

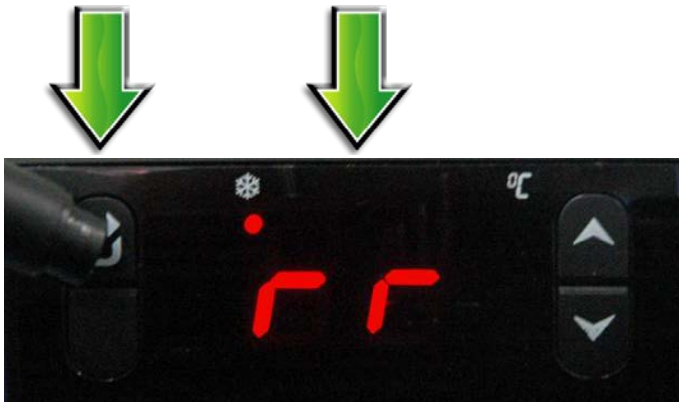

If you would like any further clarification on any of these matters please get in touch with us.

Rollex Medical NZ Limited

Postal Address: PO Box 204159, Highbrook, Auckland 2161, New Zealand

Physical Address: 38E Highbrook Drive, East Tamaki, Auckland 2013, New Zealand

Freephone: 0800 249 3722

Phone: +64 9 274 0707

Fax: +64 9 274 0709

Email (Sales): nzsales@rollexmedical.co.nz

Email (Service): nzservice@rollexmedical.co.nz