



# MATOS<sup>®</sup> WIRELESS EXTERNAL SENSOR INSTALL TIPS

Dear Customer,

The ideal location for your External MWM Sensor is approximately the middle of the chamber (in most fridges this is approx. shelf 3 or 4).

We recommend the sensor is affixed to the underside of a shelf to keep it from being knocked around by product being moved in and out of the fridge chamber. Care needs to be taken to ensure the sensor tip is not physically making contact with the metallic shelf, or the walls of the chamber.

We recommend the sensors should be held in place with something simple like a bulldog clip, as this allows for moving the sensor around the fridge without too much trouble. Normally the sensor is moved between shelves once a week. A rubber band can also easily be used for this purpose.

**Do also note the MWM Sensor should also not be touching any stock as that can give inconsistent readings.**

## Bulldog Clip Example:



## Rubber Band Example:



Please call your Rollex Medical Office if you have any questions.

### ROLLEX MEDICAL NZ LTD

PO Box 204159, Highbrook, Auckland 2161, New Zealand  
Phone: +64 9 274 0707 Fax: +64 9 274 0709  
Email: [sales@rollexmedical.co.nz](mailto:sales@rollexmedical.co.nz) [www.rollexmedical.co.nz](http://www.rollexmedical.co.nz)

### ROLLEX MEDICAL (AUST) PTY LTD

PO Box 170 | Mt Kuring-gai 2080 | Sydney, NSW | Australia  
Phone: 1300 880 441 | Fax: 1330 880 451  
Email: [sales@rollexmedical.com](mailto:sales@rollexmedical.com) | [www.rollexmedical.com.au](http://www.rollexmedical.com.au)